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Textiles are coated mostly by spreading the coating paste regularly onto the carrying material by means of a doctor knife. To also enable laboratory tests we have developed this hand-coating device which consists of the following integral parts:

1. Guiding and supporting frame
2. Coating head
3. Pin frame (material holder)

Range of Application

The coating device enables practically every coating method like:

- Floating knife
- Roller knife
- Rubber blanket knife

A number of accessories enable not only knife application, but also:

- Laminating
- Roller application

1. Guiding and Supporting Frame

The guiding shafts and bearings for the motion in longitude are accommodated in this frame.

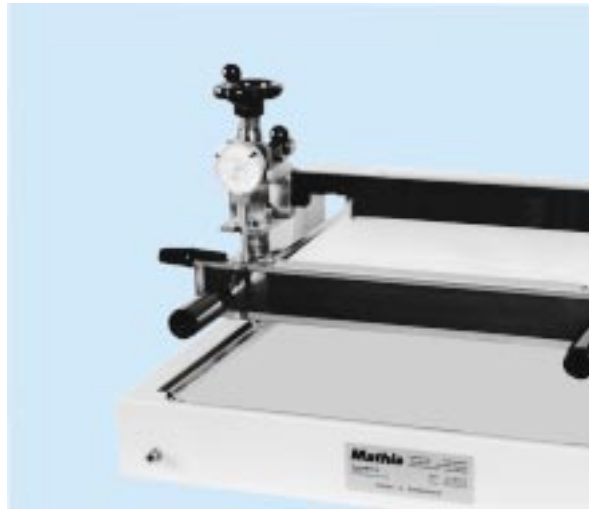
2. Coating Head

Depending on the rollers or rubber blanket employed, it is possible to work either with air knife, roller knife or rubber blanket knife.

The doctor knife is added vertically, the angle position being adjustable. The coat thickness can be adjusted and read off from measuring clocks (accuracy 0.01 mm).

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
By means of two handles the coating head is guided over the carrying material.

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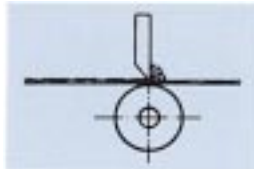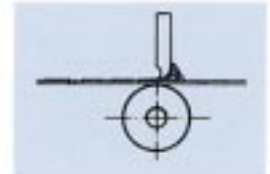
Different knife shapes

Suitable coating knives are available for PVC and PU coatings.

form B
suitable for PVC

Generally

The coating device type SV can be used in combination with our laboratory steamer/driver type DH (also DHe), since the material carrier (for paper and fabric) is suited for both machines, so that the coated material can be thermically treated without any re-pinning.

Our manufacturing programme includes:

Laboratory-Padder vertical, horizontal
Drying and fixing unit
Thermolising range
Steamer – HT-steamer
Pad-steam range
Pressure steamer
Winch
Jet dyeing apparatus
Jigger

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3. Material holders

Pin frame adjustable in length

to stretch material which can be pinned up. The tension is adjustable only in length and can be read off on a scale.

sample size: max. 33 x 43 cm

Pin frame adjustable in length and in width

to stretch material which can be pinned up. The tension is adjustable in length and in width.

The pin frames are also available with longer needles, which make it easier to hold carpet samples.

Paper and foil frame

to hold release paper, foils, non-woven fabrics etc. This can be done by means of a clamping device. The sample is tensioned in length.

Reversing pin frame

on this special pin frame the material is tensioned in length. The inner part of the frame with the pinned material can be turned. In this way it is possible to coat both sides of the sample without re-pinning.

Sieve frame

the sieve frame made of stainless steel wire netting with removable cover is specially suited to place loose material.

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4. Accessories

Laminating device type SV-K

This additional part consists of a holding device with pressing roller plus material carrier. The laminating pressure is adjustable. This device is placed at the coating head instead of the doctor blade.

Application: laminating of carriers after coating of the adhesive e. g. PVC, dispersions etc. in the reversing method.

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Fast laminating device type K-SK

Essentially the same as type SV-K however with a fast lifting device of the laminating roller. Laminating is done against a separate roller. The complete fast laminating device is to be placed behind the doctor knife and fastened by means of a clamping lever. This enables a laminating of the carrier immediately after spreading of the adhesive. Application: laminating of carriers immediately after coating of the adhesive, especially solvent containing adhesives.

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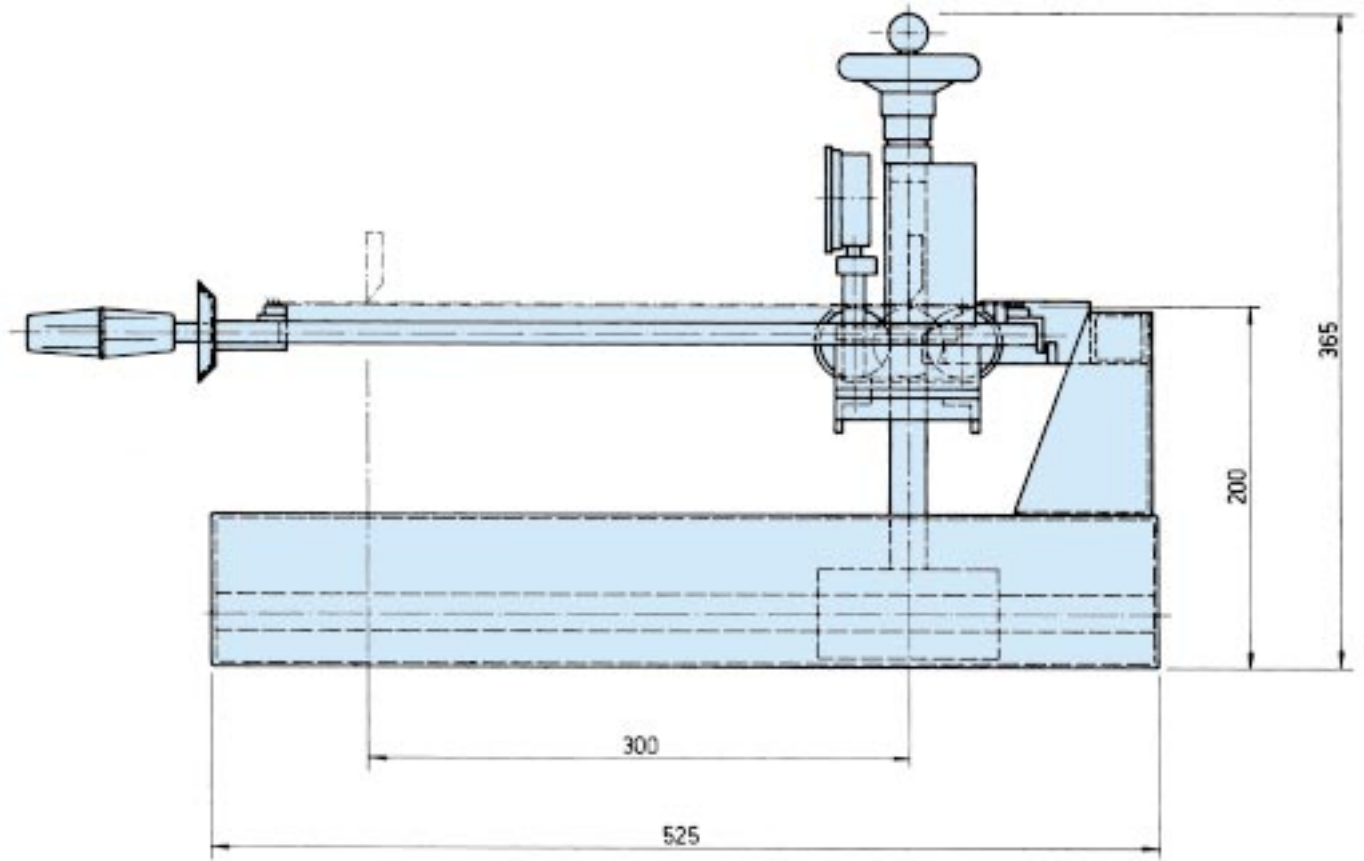
Roller coating device type SV-WB

For carpet back coatings, coatings of velvet etc. by means of roller application. The coating roller is mechanically driven. There are two doctor blades, one for the applying roller and the second one for the carpet.

The roller coating device is simple attachable. It is placed onto the holding device instead of the supporting rollers.

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max. 33 x 43 cm
 max. 28 x 30 cm
 ca. 30 kg netto
 max. 73 cm

Max. size of substrates:
 Max. size of coated sample:
 Weight:
 Width:

33 x 43 cm
 28 x 30 cm
 approx. 30 kg net
 max. 73 cm

Textilmaschinen Textile Machines
 Laborapparate Laboratory Equipment
 Beschichtungsanlagen Coating Ranges

Werner Mathis AG
 Rütisbergstrasse 3
 CH-8156 Oberhasli/Zürich
 Switzerland

Telefon 41(0)1-852 5050
 Telefax 41(0)1-850 6707
 E-mail info@mathisag.com
 Homepage www.mathisag.com

Tochterfirmen / sister companies:

Werner Mathis U.S.A. Inc.
 2260 HWY 49 N.E. / P.O. Box 1626
Concord N.C. 28026, U.S.A.
 Phone: 1-704-786-6157
 Fax: 1-704-786-6159

Aparelhos de laboratorio Mathis Ltda
 Estrada dos Estudantes 261
CEP 06700 COTIA S.P., Brasil
 Phone: 55 (0)11-7922-2333
 Fax: 55 (0)11-7922-2598
 E-mail mathbra@ibm.net